

Protocol: Target Enrichment

Exome Capture of ThruPLEX® Libraries with Illumina Nextera® Rapid Capture Exome Enrichment Kit

Required Reagents

ThruPLEX library preparation kit (one of the following required):

- ThruPLEX DNA-seq Kit (Rubicon Genomics)
 - 12 rxns, 6 single indexes (CAT. NO. R400523)
 - 12 rxns, 12 single indexes (CAT. NO. R400429)
 - 48 rxns, 12 single indexes (CAT. NO. R400428)
 - 48 rxns, 48 single indexes (CAT. NO. R400427)
 - 48 rxns, 48 dual indexes (CAT. NO. R400406)
 - 96 rxns, 96 dual indexes (CAT. NO. R400407)
- ThruPLEX Plasma-seq Kit (Rubicon Genomics)
 - 12 rxns, 12 single indexes (CAT. NO. R400490)
 - 48 rxns, 48 single indexes (CAT. NO. R400491)
 - 96 rxns, 96 dual indexes (CAT. NO. R400492)

Blocking oligos (both required):

- xGen® Universal Blocking Oligo – TS HT-i5 (Integrated DNA Technologies)
- xGen Universal Blocking Oligo – TS HT-i7 (Integrated DNA Technologies)

Illumina Capture Reagents (One of the following required):

- Nextera Rapid Capture Exome Enrichment Kit (FC-140-1000; FC-140-1001; FC-140-1002; FC-140-1003; FC-140-1083*; FC-140-1086; FC-140-1089).
*This configuration was used for protocol development
- Nextera Rapid Capture Expanded Exome Enrichment Kit

Other Consumables and Equipment

Refer to Appendix A “Consumables and Equipment” (page 44-46) in the Nextera Rapid Capture Enrichment Reference Guide, Catalog # FC-140-9001DOC, Part # 15037436 Rev. J

ThruPLEX Library Preparation

1. Prepare ThruPLEX libraries
 - Follow Section D in the ThruPLEX® DNA-seq Kit or ThruPLEX Plasma-seq Kit Instruction Manual
2. Perform Library Purification by AMPure® XP beads
 - Follow Section E.V. in the ThruPLEX DNA-seq Kit or ThruPLEX Plasma-seq Kit Instruction Manual
 - **CAUTION:** For the final elution, DNA must be eluted by re-suspending the beads in 30µL of PCR grade water, not TE Buffer

Capture of ThruPLEX Libraries

1. Combine 500 ng of each uniquely indexed ThruPLEX library
 - NOTE: The number of libraries that may be pooled is determined by the kit configuration of the Nextera Rapid Capture Exome Enrichment kit purchased (1-plex to 12-plex)
2. If necessary, adjust the volume of the pooled libraries to 38 µl
 - If the total volume of libraries is >38 µl, use a vacuum concentrator without heating to reduce volume to 38 µl
 - If the total volume is <38 µl, increase the volume to 38 µl with nuclease free water
3. Combine the following in a 0.2 ml PCR tube:
 - 38 µl DNA Library or Sample Pool
 - 1 µl xGen Universal Blocking Oligo – TS HT-i5
 - 1 µl xGen Universal Blocking Oligo – TS HT-i7
 - 50 µl buffer EHB (from Nextera kit)
 - 10 µl capture oligos CEX, EEX, or RCO (from Nextera kit)
4. Continue with Procedure step 2, page 18 of the probe hybridization per the Nextera Rapid Capture Enrichment Reference Guide, Catalog # FC-140-9001DOC, Part # 15037436 Rev. J, “Shake at 1200 rpm...”
5. For Procedure step 1 of the Second Hybridization (page 22) add the following to each well containing 25 µl of product from the first capture:
 - 13 µl buffer RSB (From Nextera kit)
 - 1 µl xGen Universal Blocking Oligo – TS HT-i5
 - 1 µl xGen Universal Blocking Oligo – TS HT-i7
 - 50 µl buffer EHB (from Nextera kit)
 - 10 µl capture oligos CEX, EEX, or RCO (from Nextera kit)
6. Continue with Procedure step 2, page 22 of the probe hybridization per the Nextera Rapid Capture Enrichment Reference Guide, Catalog # FC-140-9001DOC, Part # 15037436 Rev. J, “Shake at 1200 rpm...”

ThruPLEX® is a registered trademark of Rubicon Genomics, Inc.
Nextera® is a registered trademark of Illumina, Inc.
xGen® is a registered trademark of Integrated DNA Technologies, Inc.
AMPure® is a registered trademarks of Beckman Coulter, Inc.

ThruPLEX® Kits are intended for **Research Use Only**. It may not be used for any other purposes including, but not limited to, use in diagnostics, forensics, therapeutics, or in humans. ThruPLEX may not be transferred to third parties, resold, modified for resale or used to manufacture commercial products without prior written approval of Rubicon Genomics, Inc. ThruPLEX DNA-seq and ThruPLEX Plasma-seq are protected by U.S. Patents 7,803,550; 8,071,312; 8,399,199; 8,728,737 and corresponding foreign patents. Additional patents pending.